

MASTERflex[®]



Quick Start **Guide**

1 Go to masterflexlive.coleparmer.com or download the app from the iTunes Store or Google Play.

2



The screenshot shows the MasterflexLive login page. At the top center is the logo "MASTERFLEXLIVE™" in blue. Below it is the text "Welcome back! Please log in to your account." There are two input fields: "Email" with the text "PumpUser@coleparmer.com" and "Password" with a masked password of seven dots. A "Forgot Password?" link is located to the right of the password field. At the bottom, there are two buttons: a blue "Sign In" button and a white "Create Account" button. The "Create Account" button is highlighted with a dashed green border, and a green arrow points from a green-bordered box below it containing the text "Enter info and click **Create Account**".

MASTERFLEXLIVE™

Welcome back! Please log in to your account.

Email
PumpUser@coleparmer.com

Password
●●●●●●●

Forgot Password?

Sign In Create Account

Enter info and click **Create Account**



Register an Account

Company *	Email *
Cole-Parmer	PumpUser@coleparmer.com
First Name *	Last Name *
Test	Account 1
Country Code *	Phone Number *
United States +1	(xxx)-xxx-xxxx
Password	Password Confirmation
●●●●●●	●●●●●●

Password must contain at least 6 characters on lowercase, one numeric and one special character.



Already have an account? [Sign In](#)

Enter info and click **Register**



You have successfully created an account, please sign in with your new account.



Back to Sign In



Click **Back to Sign In**



Welcome back! Please log in to your account.

Email

PumpUser@coleparmer.com


Password

●●●●●●●●

[Forgot Password?](#)



Enter your info and click **Sign In**

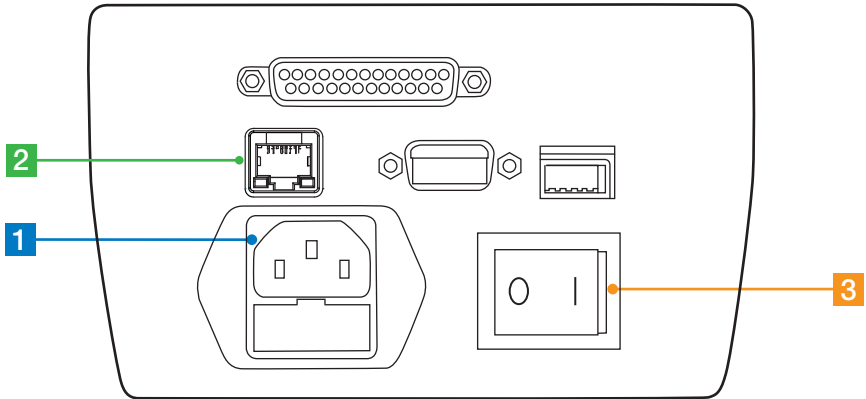

*Diagram for illustration only, your unit may be different

Step 1

Plug in power cord

Step 2


Connect drive to an internet-ready network using the RJ-45 connection port on the rear of your device or with wifi enabled capability.

Step 3


Power on the drive- drive is now ready to connect to your account.

Page Navigation

Click to view/hide
the left hand
menu bar

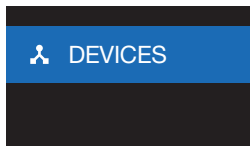

Select **Account Settings**,
My Profile or **Sign Out**
from the drop down menu

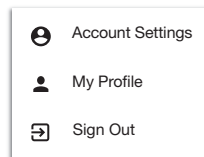

For more information on how to add a device to your MasterflexLive account and how to access the pump controls, please visit masterflex.com/masterflexlive.

MASTERFLEX[®]

US & Canada only

Toll Free 1-800-MASTERFLEX | 1-800-637-3739

Outside US & Canada

***EN809 manufactured by:**

Cole-Parmer Instrument Company

28W092 Commercial Avenue, Barrington, IL 60010

A-1299-5203 Edition 01